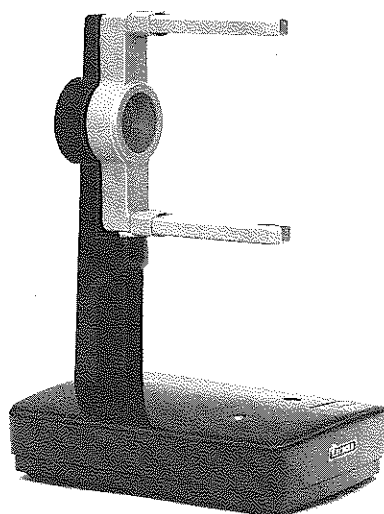
Experimentierständer Helmholtz-Spulenpaar

Universal Stand Pair of Helmholtz-Coils

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555 06

Fig. 1

The universal stand has been designed for mounting the electronic tubes (555 07 to 555 17) and the optical analogue (555 19).

The pair of Helmholtz coils is fixed on the universal stand. It is used for experiments with the electronic tubes (555 07, 555 10 to 555 12 and 17). The coils produce a homogenous magnetic field of variable strength, in which the deviation of electrons is shown and analysed.

Notes

1. The five-digit numbers quoted in brackets refer to the catalogue numbers of the respective apparatus.
2. The specifications and illustrations are not binding in every detail for the design of the apparatus. It is our policy always to keep our manufacturing programme right up to date so that it makes full allowance for the developments acquired in all scientific and technical fields.

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A. Universal Stand (555 05)

1. Description

The large base-plate (1) (dimensions: approx. $240 \times 180 \times 50 \text{ mm}^3$) and the high stanchion (2) (approx. 280 mm high), are made of a light metal alloy, hot-enamelled in blue. The handle is covered with black imitation leather. Three rubber feet assure a non-skid mounting.

Two arms (3) (about 160 mm long and spaced 140 mm apart) hold the various tubes. The arms are made of a plastic material, which has excellent temperature and electrical insulation features. These arms, i. e. the hoop, are provided with two recesses to fix the tubes at their plastics caps. The tubes can be locked in place by means of two sliders (4). The arms are mounted on the stanchion and can be rotated about a horizontal axis so that the tubes can be operated in any desired position about their longitudinal axis (Fig. 2).

The two holes (5) in the base-plate serve to fix the Helmholtz coils. In the stanchion there is a 4 mm dia earthing socket (6).

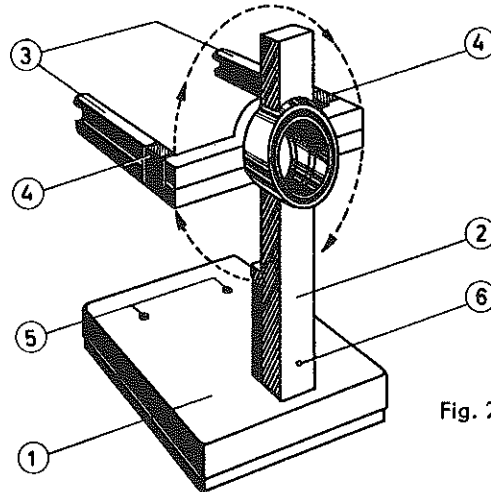
Fig. 2

2. Use

The arms (3) are turned in the required position.

To fix the tubes at both ends, they are pushed through the opening of the stanchion. When the sliders (4) are moved back, the plastics caps are fitted in the recesses of the arms. It is recommended to spread the arm with one hand, while the other hand inserts carefully the plastics caps of the tube into the provided recesses. **In no case** the cap must be used to force the arm apart! When releasing the tube, the same method has to be applied.

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After having fixed the tube, it is locked in place by means of the sliders.

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1. Description

Each of the coils is wired on a plastics coil-former of 136 mm average \varnothing . Each coil has 320 windings. The coil-former is mounted on a holding rod, (approx. 140 mm long) on the other end of which there is an insulating button. Two rifflled sliding plastics sleeves serve to mount the holding rods on the universal stand in an insulating manner.

Both ends of the wire-winding are connected to the 4 mm plug, which is laterally fixed to the coil-former or to the holding rod. Within the holding rod there is a 4 mm connection socket.

The operating current of the coils is 1 to 1.5 A, corresponding to an operating voltage of approx. 10 to 15 V, if connected in series.

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2. Use

After having clamped the tube in question in vertical position of the arms, the coils are inserted with their holding rods into the holes of the baseplate (Fig. 3). The insulating button must fit into the provided recess. Now the holding rod is fixed by pressing the plastics cap into the tapering hole.

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$$a = 3,4 \text{ cm}$$

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$$H = 33,8 \cdot 10^2 \text{ m}^{-1} \cdot I$$

The direct current part of the three-phase current supply (522 48), the rectifier block (522 46) which is operated at the three-phase current mains, or two nickel-cadmium accumulators (522 71), each of which provided with a rheostat of 11Ω (537 26), are a suitable current source for the magnetization.

The connection is executed according to the usual series-connection of the Helmholtz-mounting via the plugs of the coil-formers. The sockets within the holding rods are connected to each other. Now the current flows through both coils in the same sense.

3. Calculation of the magnetic field

From the Biot-Savart relation it follows for the field strength H of the (nearly) homogeneous magnetic field of the pair of Helmholtz coils:

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